

6. Start BV Workbench and Set up software license

1. Click [Brainvision]-[BV Workbench] registered in the start menu or double-click the shortcut created on the desktop to start BV Workbench.



Icon of BV Workbench

***BV Workbench can be started even when MiCAM03 power is off.
In this case, it will be in the offline mode where you can load saved data and analyze data.***

2. BV Workbench starts.



3. After installing BV Workbench, the following screen will be displayed the first time you start it. After inputting/selecting information required for license issuance on this screen, click the "SEND" button to start email software.

Check information displayed in the email and send it to us so that we can receive the information. Based on that, we will create a license and we will send you the license file by email.

Please fill out the form below to request a new license.

License status : License file not found.

Name : Taro Yamada

Organization : Brainvision Inc.

E-mail : yamada.taro@brainvision.co.jp

License type : Standard

Hardware ID : BCBB-DC95-E2A1-67E1-ADEC-8DCC-927F-AAF7/5161-B295-5D1F-4B6B-5F7D-FF72-E24B-2061/4490-271C-3A5B-5266-BC7F-4CB7-DB1C-A378/26BF-A695-9E6F-4AD4-9561-C4E3-984F-AAF5/FBD5-E830-5F5E-91C3-A9BA-1214-63D6-8A37/ECC1-A393-8FDC-EA3E-29DE-C5AB-5C2A-AABF

Required features

- Acquisition (to disk) : Acquire data directly to disk (for very long acquisitions)
- Acquisition (time lapse) : Acquire data in time lapse mode
- ESTM-8 : For ESTM-8 users
- ESTM-9 : For ESTM-9 users
- MiCAM03 : For MiCAM03 users
- MiCAM05 : For MiCAM05 users
- Third party camera : For third party camera users
- Post processing : Image processing module (filters)
- Analysis : Analysis module
- E-Field : Electric field imaging module

Click to start email software

BV Workbench licenses are issued for each computer by using a hardware ID created from information unique to an installed computer.


Licenses issued for other computers cannot be used.

Please contact SciMedia Ltd. (inquiry@scimedia) for reissuance of license.


4. Set up a license in BV Workbench using a license file sent from us.

Again, click [Brainvision]-[BV Workbench] registered in the start menu or double-click the shortcut created on the desktop to start BV Workbench.


The following window will open. Click [CLOSE] button.

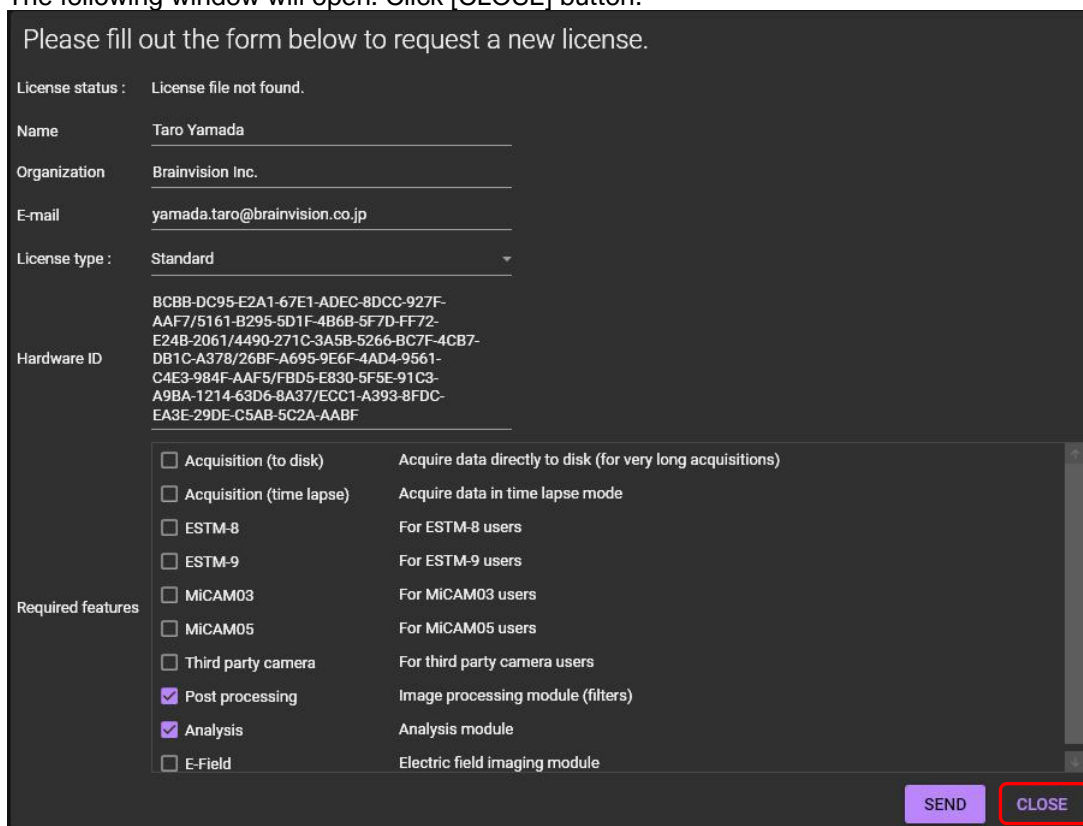

5. Click the following icon (“App settings”) on the toolbar to open the [App settings] screen.

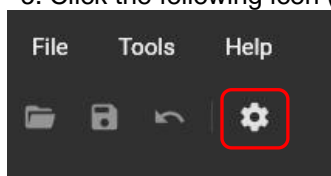

6. Click [IMPORT LICENSE FILE...] at the bottom of the [App settings] screen to display the file selection screen.

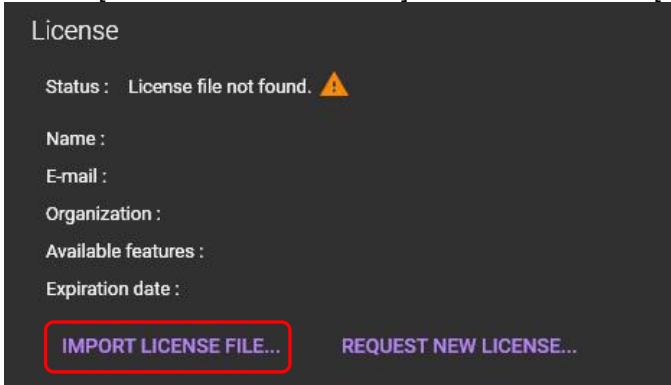

7. On the file selection screen, select a license file (extension: “lic”) sent by us and click the [Open] button.
When the license is set, the license information is displayed as shown below and this software can be used.

